



Dricyclers



Applications

- Storing reagents or soil samples
- Drying biological specimens
- Cooling crucibles
- Storage sludge samples

Features

- Durable, hand crafted construction. Fabricated from stainless steel and glass
- Superior seal integrity. One piece neoprene door gasket
- Rigid door with full length door hinges
- Shelves are adjustable in half-inch increments
- Two different sizes to select from
- Automatic desiccant regeneration available
- Automatically regenerates the silica gel desiccant enclosed in the control unit eliminating the need to replace or bake desiccant
- Easy to read door mounted hygrometer.

Boekel Dricyclers Drying Sequence

Dricyclers automatically regenerate desiccant every five hours.

4.5 hours - Silica gel cassette housed in control box is exposed to desiccator air for 4.5 hours.

0.5 hour - Silica gel cassette closed off from desiccator air and regenerated by heating. Moisture is vented to outside air.

Specifications

Model	Model 134041 (-2)	Model 134241 (-2)	Model 134441 (-2)
Overall W x L x H	40.6 cm x 31.4 cm x 31.1 cm (16" x 12 1/2" x 12 1/4")	40.6 cm x 31.4 cm x 31.1 cm (16" x 12 1/2" x 12 1/4")	52.1 cm x 58.4 cm x 57.2 cm (20 1/2" x 23" x 22 1/4")
Chamber W x L x H	29.2 cm x 25.4 cm x 30.5 cm (11 1/2" x 10" x 12")	29.2 cm x 25.4 cm x 30.5 cm (11 1/2" x 10" x 12")	40 cm x 50.2 cm x 55.2 cm (15 3/4" x 19 3/4" x 21 3/4")
Net Weight	11.8 kg (26 lbs)	11.8 kg (26 lbs)	27.2 kg (60 lbs)
Power	115VAC (or 230V) 50/60 Hz 35 W	115VAC (or 230V) 50/60 Hz 35 W	115VAC (or 230V) 50/60 Hz 35 W

Ordering Information

Description	Catalog/Model Number 115V	Catalog/Model Number 230V
Dricycler, small, 2 ceramic shelves,	134041	134041-2
Dricycler, small, 2 stainless steel shelves	134241	134241-2
Dricycler, large, 2 stainless steel shelves	134441	134441-2

Accessories

Shelves' Catalog Number	Use with
C000227	134041, 134041-2
D1004369	134241, 134241-2
D1004371	134441, 134441-2