

Microscope Slide Cabinet Dolly Safety Cautions

- Never sit on, stand on, or otherwise ride the dolly.
- Before loading/unloading the dolly, ensure all casters are locked.
- Follow ergonomically sound practices when loading/unloading the dolly.
- Never stack the dolly higher than 10 microscope slide cabinets (< 5ft or 1.5m tall)
- If the dolly is used without a cabinet base, ensure bottom cabinet is properly latched to dolly.
- Load heaviest (fullest) microscope slide cabinets on the bottom.
- Ensure all stacked cabinets are properly connected to each other before moving the dolly.
- Load cabinets drawers from the bottom up. Never load slides in the top cabinets first.
- Do not open more than one cabinet drawer at a time, as this increases the risk of toppling.
- Push loaded dolly from the center, never from the top.
- Avoid walking backwards/pulling the loaded dolly.
- Lock all casters when the dolly is not in use (i.e. dolly/cabinet tower is in storage location).
- Do not use for other purposes than intended.
- Do not use on staircases or on ramps.
- Do not use in any manner that may result in personal injury or damage to property.



**TIP OVER
HAZARD**



**Warning
Toppling
hazard**