

PRODUCT INFORMATION

Boekel's line of Stainless Steel & Glass Desiccators are designed as premium quality, air tight, moisture proof cabinets. The desiccators provide excellent storage environments for materials and parts requiring storage at low humidity. The glass and stainless steel combination provides a sturdy stackable cabinet which allows the user to see what is being stored without opening the door and exposing the material to room humidity. The cabinet door is fully gasketed and comes with a very secure latching mechanism.

Each Boekel desiccator comes with two shelves and a single tray for holding desiccant. The model 1340 has two 9 3/4" x 8" stainless steel shelves, the model 1344 has two 19" x 13 1/2" stainless steel shelves. Shelf separation on the model 1344 is a minimum of 3/4". All models come with a manual relief valve installed on the upper right side of the unit.

Optional additional shelving and replacement door gaskets are available for all models.

CAUTION

Boekel's Desiccators are not intended for the purpose of maintaining a vacuum or pressure.

UNPACKING

1. Remove desiccator from carton.
2. Inspect for damage. Report all shipping damage to the carrier immediately. Shipping damages are covered by the carrier and repair and/or replacement for shipping damage must be coordinated through the carrier.
3. Retain all packaging material in the event that the unit must be returned.
4. Completely fill out and return the self addressed Boekel Warranty Registration card to Boekel Scientific.

SETUP

Place the desiccator on a flat surface and install the two shelves. Desiccator models 1340 and 1342 may be stacked up to four units high without external bracing. Desiccator model 1344 may be stacked two units high without bracing. Stacking may be limited by the total weight of the objects to be stored in the desiccator.

CLEANING

1. The unit interior and exterior may be cleaned with a damp cloth or sponge and warm water.
2. If an internal spill occurs, swab interior of unit with a suitable bactericide or disinfectant.
3. When cleaning door, use a soft cloth and water or mild detergent.

NOTE: Abrasive materials such as paper towels may scratch door gasket. **DO NOT USE SOLVENTS.**

SERVICE: RETURN AUTHORIZATION REQUIRED.

Should service be required, contact your salesperson or call Boekel Scientific at 215-396-8200 or 800-336-6929.

SPECIFICATIONS

SMALL DESICCATOR

Models 1340 and 1342

Exterior: 12 1/2"W x 12 3/8"H x 12 3/8"D

LARGE DESICCATOR Model 1344

Exterior: 16 3/4"W x 22 1/4"H x 23"D

Maximum Temperature: 150°C