# **OPERATING INSTRUCTIONS**



## DRICYCLER DESICCATORS

MODELS 134041, 134041-2, 134241, 134241-2, 134441 & 134441-2

### PRODUCT INFORMATION

Boekel's Dricycler Desiccators are designed to eliminate the need to manually change desiccant on a periodic basis. The Dricycler will establish and maintain on average, a relative humidity in the 25% to 35% RH range without the addition of desiccants. CAUTION: These Dricycler Desiccators are not intended for the purpose of maintaining a vacuum or pressure.

#### **UNPACKING**

- 1. Remove the Dricycler from carton.
- 2. Check to see that the Dricycler Desiccator was not damaged during shipment.
- 3. Retain all packing material in the event that the unit must be returned.
- 4. Completely fill out and return the self-addressed Boekel warranty card.

#### **ASSEMBLY**

There is no assembly required for the Dricycler Desiccator. Simply remove it from the box and plug it in to an appropriate outlet.

## **OPERATING INSTRUCTIONS**

- Plug unit into the appropriate power outlet
- 2. The green Power indicator light should immediately go on.

SPECIFICATIONS: SMALL DRICYCLERS

DIMENSIONS: 16 1/2" X 12 3/8" X 12 3/8"

POWER: 115 VAC (35 W) 134041

115 VAC (35 W) 134241 230 VAC (35 W) 134041-2 230 VAC (35 W) 134241-2

Maximum Temperature: 150°C

- 3. The red indicator will occasionally go on indicating that the regenerating auto dry process is active.
- 4. Upon initial start-up, let the Dricycler run for a period of 24 hours before loading. This allows stabilization of the unit chamber.

#### **CLEANING**

- Disengage power cord before performing any cleaning.
- 2. The unit interior and exterior may be cleaned with a damp cloth or sponge and warm water.
- If an internal spill occurs swab interior of unit with a suitable bactericide or disinfectant.
- 4. When cleaning the door, use a soft cloth and water or mild detergent.

NOTE: Abrasive material such as paper towels may scratch the door or gasket. DO NOT USE SOLVENTS,

SERVICE: RETURN AUTHORIZATION IS REQUIRED.

Should service be required, contact your salesperson or call Boekel Scientific at 215-396-8200 for return authorization.

SPECIFICATIONS: LARGE DRICYCLERS

DIMENSIONS: 20 1/2" X 22 1/4" X 23" POWER: 115 VAC (35 W) 134441 230 VAC (35 W) 134441-2

Maximum Temperature: 150°C