

DRICYCLER DESICCATORS

MODELS 134041, 134041-2, 134241, 134241-2, 134441 & 134441-2

PRODUCT INFORMATION

Boekel's Dricycler Desiccators are designed to eliminate the need to manually change desiccant on a periodic basis. The Dricycler will establish and maintain on average, a relative humidity in the 20% to 35% (dependent upon the ambient room humidity). Without the addition of desiccants. **CAUTION:** These Dricycler Desiccators are not intended for the purpose of maintaining a vacuum or pressure.

UNPACKING

1. Remove the Dricycler from carton.
2. Check to see that the Dricycler Desiccator was not damaged during shipment.
3. Retain all packing material in the event that the unit must be returned.
4. Completely fill out and return the self-addressed Boekel warranty card.

ASSEMBLY

There is no assembly required for the Dricycler Desiccator. Simply remove it from the box and plug it in to an appropriate outlet.

OPERATING INSTRUCTIONS

1. Plug unit into the appropriate power outlet and turn power switch to "ON" position.

SPECIFICATIONS: SMALL DRICYCLERS

DIMENSIONS: 16 1/2" X 12 3/8" X 12 3/8"

POWER: 115 VAC (35 W) 134041
115 VAC (35 W) 134241
230 VAC (35 W) 134041-2
230 VAC (35 W) 134241-2

Maximum Temperature: 150°C

2. Upon initial start-up, let the Dricycler run for a period of 24 hours before loading. This allows stabilization of the unit chamber.

CLEANING

1. Turn power switch to "OFF"
2. Disengage power cord before performing any cleaning.
3. The unit interior and exterior may be cleaned with a damp cloth or sponge and warm water.
4. If an internal spill occurs swab interior of unit with a suitable bactericide or disinfectant.
5. When cleaning the door, use a soft cloth and water or mild detergent.

NOTE: Abrasive material such as paper towels may scratch the door or gasket. **DO NOT USE SOLVENTS,**

SERVICE: RETURN AUTHORIZATION IS REQUIRED.

Should service be required, contact your salesperson, or call Boekel Scientific at 215-396-8200 for return authorization.

SPECIFICATIONS: LARGE DRICYCLERS

DIMENSIONS: 20 1/2" X 22 1/4" X 23"

POWER: 115 VAC (35 W) 134441
230 VAC (35 W) 134441-2

Maximum Temperature: 150°C